

# Quick Guide


## Poultry



### Step 1

#### Insert the sample chamber

Insert the sample chamber in the direction indicated by the arrow.

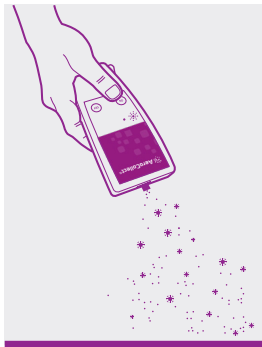

### Step 2

#### Start! Press ON (green)


The instrument can be started by pushing the green ON-button.



### Step 3

#### Kick up dust from the floor


Kick up the bedding material to facilitate more airborne particles.

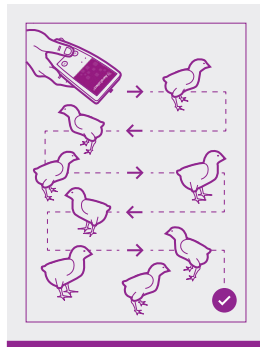

### Step 4

#### Sampling from entire house


Walk a route in the entire house for a minimum of 5 minutes to ensure a reliable result.

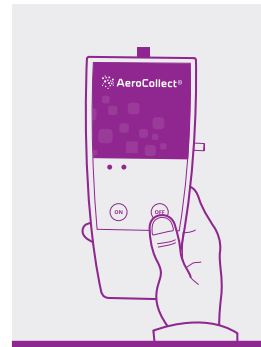

### Step 5

#### Stop! Press OFF (red)

The sampling procedure is stopped by pressing the red OFF-button.

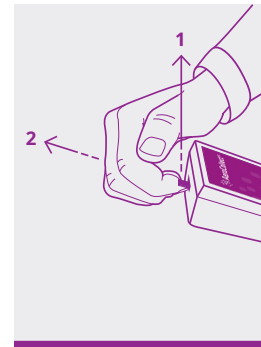

### Step 6

#### Remove the sample chamber

Lift up (1) the front of the sample chamber while pulling it out (2) of the AeroCollect®.





Watch how to use AeroCollect®  
Scan the QR code with the camera on your phone



# Quick Guide

## Poultry

